



FIREPol[®] DNA Polymerase

Cat. No.	Pack Size	Conc.
01-01-0000S	100 U SAMPLE	5 U/ μ l
01-01-00500	500 U	5 U/ μ l
01-01-01000	1000 U	5 U/ μ l
01-01-02000	2000 U	5 U/ μ l

For *in vitro* use only

Description:

FIREPol[®] is a highly processive, thermostable DNA polymerase. Due to its genetic modifications FIREPol[®] has an enhanced stability at room temperature with no activity loss for up to 1 month. The enzyme has 5'→3' polymerase and 5' flap endonuclease activity.

Source:

Purified from an *E. coli* strain that carries an overproducing plasmid containing a modified gene of *Thermus aquaticus* DNA Polymerase.

Applications:

- Suited for a wide range of PCR assays
- TA cloning

Reagents Provided:

- **FIREPol[®] DNA Polymerase**
- **10x Reaction buffer B** (Mg²⁺ free)
0.8 M Tris-HCl, 0.2 M (NH₄)₂SO₄, 0.2% w/v Tween-20
- **10x Reaction buffer BD** (Mg²⁺ and detergent free)
0.8 M Tris-HCl, 0.2 M (NH₄)₂SO₄
- **25 mM MgCl₂**
- **10x Solution S**
Additive that facilitates amplification of difficult templates (e.g. GC-rich DNA templates). This solution should be used at a defined working concentration (1x, 2x or 3x solution).
Solution S is NOT a reaction buffer and should be used ONLY IF non-specific amplifications occur.

Concentration:

5 U/ μ l

Unit definition:

One unit is defined as the amount of enzyme required to catalyze the incorporation of 10 nmol of dNTPs into an acid-insoluble form in 30 minutes at 74°C.

Storage and Dilution buffer:

50% glycerol (v/v), 20 mM Tris-HCl pH 8.7 at 25°C, 100 mM KCl, 0.1 mM EDTA and stabilizers.

Quality control:

The enzyme is free of nicking and priming activities, exonucleases and non-specific endonucleases. SDS/PAGE - 95 kD band, >98% pure. Activity and stability tested via thermo-cycling. The error rate per nucleotide per cycle is ~ 2.5 x 10⁻⁵; the accuracy is ~ 4 x 10⁴. Estimated half life at 95°C is 1.5 hours.

Shipping and Storage conditions:

Routine storage: -20°C

Shipping and temporary storage for up to 1 month at room temperature has no detrimental effects on the quality of FIREPol[®] DNA Polymerase.

Recommended PCR reaction mix:

Component	Volume	Final conc.
FIREPol [®] DNA Polymerase (5 U/ μ l)	0.4-1.0 μ l	0.02-0.05 U/ μ l (2-5 U)
10x Buffer B or BD	10 μ l	1x
25 mM MgCl ₂	6-10 μ l	1.5-2.5 mM
dNTP mix (20 mM of each)	1 μ l	200 μ M of each
Primer Forward (10 pmol/ μ l)	1-3 μ l	0.1-0.3 μ M
Primer Reverse (10 pmol/ μ l)	1-3 μ l	0.1-0.3 μ M
DNA template	variable ¹	variable ¹
10x Solution S Not for standard PCR	0, 10, 20 or 30 μ l	1x, 2x or 3x
H ₂ O PCR grade	Up to 100 μ l	
Total	100 μl	

¹Conc. of cDNA 0.01 pg/ μ l -0.1 ng/ μ l ; gDNA 0.1 ng/ μ l – 10 ng/ μ l

Recommended PCR cycles:

Cycle step	Temp.	Time	Cycles
Initial denaturation	95°C	3-5 min	1
Denaturation	95°C	15-30 s	26-35
Annealing	50-68°C	30-60 s	
Elongation	72°C	1-4 min	
Final elongation	72°C	5-10 min	1

IMPORTANT: Annealing temperature should be 2-6°C lower than the primer melting temperature. Elongation time should be ~1 min/1 kb.

In order to prevent contamination, we recommend you to setup the reaction under laminar or in PCR box.

Safety warnings and precautions:

This product and its components should be handled only by persons trained in laboratory techniques. It is advisable to wear suitable protective clothing, such as laboratory overalls, gloves and safety glasses. Care should be taken to avoid contact with skin or eyes. In case of contact with skin or eyes, wash immediately with water.

Related products:

Product name	Pack size	Cat. No.
HOT FIREPol[®] DNA Polymerase	500 U	01-02-00500
HOT FIREPol[®] DNA Polymerase	1000 U	01-02-01000
5x FIREPol[®] Master Mix (1.5 mM MgCl ₂ final conc.)	250 reactions	04-11-00115
5x FIREPol[®] Master Mix (2.5 mM MgCl ₂ final conc.)	250 reactions	04-11-00125
5x FIREPol[®] Master Mix Ready to Load (1.5 mM MgCl ₂ final conc.)	250 reactions	04-12-00115
5x FIREPol[®] Master Mix Ready to Load (2.5 mM MgCl ₂ final conc.)	250 reactions	04-12-00125
dNTP MIX (20 mM of each)	20 µmol	02-31-00020
dNTP MIX (20 mM of each)	100 µmol	02-31-00100
dNTP SET (100 mM)	4 x 25 µmol	02-21-00100
dNTP SET (100 mM)	4 x 100 µmol	02-21-00400

Permitted Use: This product is supplied for research use only (the **Permitted Use**). If the customer wishes to use the product for any purpose other than the Permitted Use, including (without limitation) resale or alteration, the customer should obtain the appropriate licence from Solis BioDyne. Some applications of this product may require a license/licenses from one or more third parties which are not provided by the purchase of this product. Users should obtain the licence if required. [Patent Notice - EP2501716 and other patents pending.](#)

Warranty and Disclaimer: This product shall comply with its relevant specification and be fit for its stated purpose, but Solis BioDyne gives no other warranty and makes no representation as to description or quality. Any such warranty or representation is excluded, to the fullest extent permitted by law. In particular, but without limiting the foregoing, Solis BioDyne shall not be liable for the failure of the product to comply with its relevant specification where such failure arises as a result of: (i) customer negligence or because the customer failed to follow any of the applicable technical data or safety sheets, standard user materials, use guidelines or any other information provided by Solis BioDyne as to the storage, transportation, handling, use or maintenance of the products or other good practice regarding the same, or (ii) the customer altering the products in any way without the prior written consent from Solis BioDyne, or (iii) the products differing from the relevant specification as a result of changes made to ensure their compliance with applicable statutory or regulatory requirements.

Nothing shall limit or exclude Solis BioDyne's liability for death or personal injury caused by its negligence, fraud or fraudulent misrepresentation or any matter in respect of which it would be unlawful for Solis BioDyne to exclude or restrict liability. Without limiting the foregoing, Solis BioDyne shall under no circumstances whatever be liable to the customer, whether in contract, tort (including negligence), breach of statutory duty, or otherwise, for any loss of profit, or any indirect or consequential loss arising under or in connection with the products and Solis BioDyne's total liability to the customer in respect of all other losses arising under or in connection with the product, whether in contract, tort (including negligence), breach of statutory duty, or otherwise, shall in no circumstances exceed the price of the products supplied in respect of which the liability has arisen.